

# Product manual

Labinco L46



LABINCO B.V.

P.O.BOX 3259-4800DG Breda, The Netherlands Zinkstraat 64-4823AC Breda, The Netherlands Tel.:0031(0)76-5417582, Fax:0031(0)76-5420173 E-mail:info@labinco-bv.com - Internet:www.labinco-bv.com



## Model L46



#### **Description:**

The model" I 46" is a vortex-mixer with strong vibrating action, for test tubes 7 mm diam. up to 35 mm. Both; "Press-to-Mix" action and "continuous" action are provided.

Stepless speed control, up to maximum speed 2500 r.p.m.

Strong and steady unit.

Lined out for accessory adaption(see <a href="http://www.labinco-bv.com">http://www.labinco-bv.com</a>).

Supplied with line-cord and standard cup.

### Operating instructions:

After carefully unpacking the vortex mixer, place the mixer on a flat table. Connect the line-cord to sufficient voltage (see specifications).

"Speed control" to be checked in ultimate clock-wise direction.

Switch" modeswitch "to" continuous" mode.

The mixer is switched "on' 'by turning "speed control' 'clock-wise, the pilot-lamp will lit and the mixer starts running. Turning "speed control" clock-wise more or less will vary mixing speed.

You can switch to "press-to-mix" mode and start mixing by pushing testtube into the rubber cup.

Specifications: 230 Volt 35 Watt 50/60 HZ

(other lines on request)

Dimensions: 135x120x135 mm (LxWxH)

Weight 3,2 KG



#### LABINCO B.V.

P.O.BOX 3259-4800 DG Breda, The Netherlands Zinkstraat 64-4823AC Breda, The Netherlands Tel.:0031(0)76-5417582, Fax:0031(0)76-5420173 E-mail:info@labinco-bv.com – Internet:www.labinco-bv.com